

VERTICAL DEEP FREEZER “MAC” MSW-135

Features

- No frost hassle free cooling system
- Microprocessor controlled digital display for temperature
- Control range -16°C to -24°C.
- Efficient and homogenous backside freezing system
- Pull out drawers
- Different size drawers for different applications
- CFC free refrigerant

Precautions

1. Use this appliance only for its intended purpose as described in this use and care guide.
2. This freezer must be properly installed in accordance with the installation instructions before it is used. See grounding instructions in the installation section.
3. Never unplug your freezer by pulling on the power cord. Always grasp the plug firmly and pull straight out from the outlet.
4. Repair or replace immediately, all electric service cords that have become frayed or otherwise damaged. Do not use a cord that shows cracks or abrasion damage along its length, the plug, or the connector end.
5. Unplug your freezer before cleaning or before making any repairs. Note: If for any reason this product requires service, we strongly recommend that a certified technician performs the service.
6. Do not use any electrical device or any sharp instrument in defrosting your freezer.
7. If your old freezer is not being used, we recommend that you remove the doors. This will reduce the possibility of danger to children.



8. After your freezer is in operation, do not touch the cold surfaces in the freezer compartment, particularly when hands are damp or wet. Skin may adhere to these extremely cold surfaces.
9. Do not leave food in this scientific deep freezer.
10. Do not use metal or any sharp object for removing the freeze ice as it can damage to surface of the freezer.

Operating Procedure

A. Opening of the door:

1. Remove the padlock if you have one.
2. Move the key into the open position by rotating it counterclockwise.
3. Grasp the latch handle (preferably with your left hand) and pull it toward yourself until the latch disengages from the cabinet strike.
4. Keep pulling by the latch handle or the integrated door handle to open the main door.

B. Closing of the door:

1. Grasp the latch handle (preferably with your left hand) and pull it toward yourself, rotating the latch into the open position.
2. Move the freezer door into the closed position and gently push the handle away from you, making sure that the latch engages fully with the cabinet strike.
3. Keep applying gentle pressure to the latch handle until the latch is securely in closed position.

C. Make sure that door is closed properly.

D. Adjust the temperature by using the arrow button.

E. Set the alarm, as per needed by using the alarm button.